

Replacement LAQUAtwin Sensor Instruction Manual


Compatibility table between sensors and meters

Measured item	Replacement sensor		Meter	
	Display on body	Model	Model	Series
pH	LAQUAtwin pH	S010	B-711, B-712, B-713	LAQUAtwin
			B-211, B-212, B-213	twin
Salt	LAQUAtwin Salt	S021	B-721	LAQUAtwin
Na ⁺	LAQUAtwin Na ⁺	S022	B-722	LAQUAtwin
K ⁺	LAQUAtwin K ⁺	S030	B-731	LAQUAtwin
NO ₃ ⁻	LAQUAtwin NO ₃ ⁻	S040	B-741, B-742, B-743	LAQUAtwin
			B-341, B-342, B-343	twin
Ca ²⁺	LAQUAtwin Ca ²⁺	S050	B-751	LAQUAtwin
Conductivity	LAQUAtwin COND	S070	B-771	LAQUAtwin

* LAQUAtwin sensors do not support CARDY series and B-173 (twin compact conductivity meter).


Replacement procedures

1. Lift the sensor tongue tip and slide the sensor a little away from the meter.
2. Pull out the sensor all the way from the meter.
3. Confirm that the waterproofing gasket is clean and undamaged.
4. Slide the sensor onto the meter so that catch "A" on the back of the meter fits into hole "a" on the sensor tongue as shown.

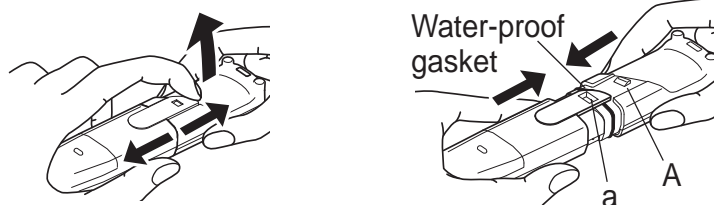

Note

- Be careful not to twist the water-proof gasket.
- Make sure that the sensor tongue is outside the meter case. If the tongue is inserted between the case and the connector of the meter, it may damage the connector.
- When removing the sensor, do not let any water get inside the meter. If some moisture remains on the waterproof gasket, wipe off it carefully.
- Perform calibration before measurement. Refer to the instruction manual of the meter for the details.

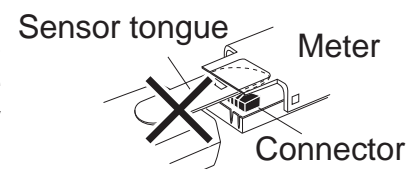

For any questions regarding this product, please contact your local agency, or inquire from the HORIBA website. (http://global.horiba.com/contact_e/index.htm)

HORIBA, Ltd.

2 Miyanohigashi, Kisshoin Minami-ku, Kyoto 601-8510 Japan

<http://www.horiba.com>